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Multi-Purpose and Analytical Balances

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Unpacking

1. Remove adhesive tape round the plastic casing.
2. Carefully unpack the balance and accessories from the two plastic halves (Fig. 1).
3. Swing aside the two bolts (A) on the weighing compartment cover and remove housing (B) (Fig. 2).
4. Unscrew bolt (C) with screwdriver.
5. Remove the two screws (E) in the weighing compartment cover and withdraw wedge (D) forwards.
6. Cut the plastic thread at the hanger and remove plastic piece (F).
7. Slacken the yellow bracket (G) with screwdriver and remove it.
8. Unscrew yellow plate (H) in housing (B) and carefully take out the beam.

Important: the two knurled nuts on the beam must not be altered.

All the packing material is now removed from the balance and setting-up can be begun.

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Setting up

1. Install the balance at the place where it is to be used.
2. Thoroughly clean the balance and accessories with a cleaning cloth. The sapphire pins, knife-edges and bearings must be scrupulously clean (see the Operating Instructions under "Maintenance").
3. Check the voltage at terminal strip ① (to change the voltage, see the Operating Instructions).
4. Arrest the balance with locking lever ⑫ (put lever ⑫ in position I) (Fig. 4).
5. Unclip damping pot cover ③ and pull downwards.
6. Carefully insert beam (Fig. 3) and replace damping pot cover ③.
7. Insert hanger plate ② and fit hanger.
8. Check that all weights ⑥ are in the grooves.
9. Remove locking lever ⑫.
10. Fit housing ⑦ and secure with bolt ⑨.
11. Fit operating knobs ⑧ and locking lever ⑫.
12. Fit pan brake ⑪ (Fig. 5).
13. Hang weighing pan on upper hook ⑤.
14. Adjust the foot screws until spirit level ⑩ reads correctly.
15. With type H 15 or 16: insert the magnifier in the front frame from left to right.
16. Insert the mains plug.

For checking and weighing see the Operating Instructions. After setting up wait for about 30 minutes until the balance has attained the ambient temperature.

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