COMMANDE D'ALIMENTATION DU LASER SPECTRA-PHYSICS STABILITE 2017

FICHE N° 5265

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Fabricant : Spectra Physics ; Spectra Physics

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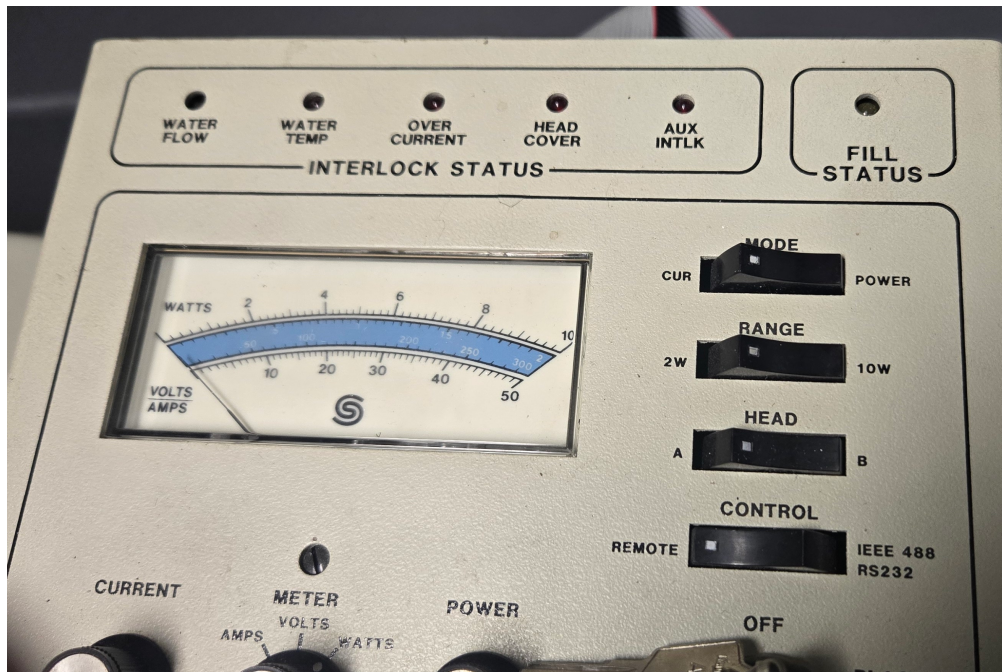
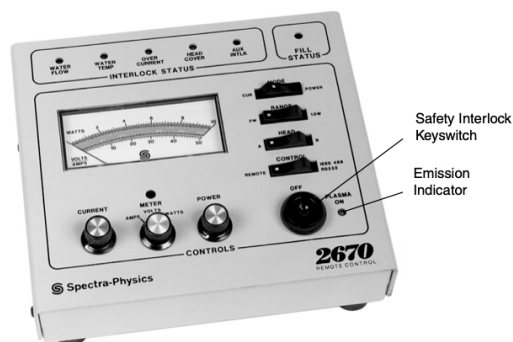
Figure 2-4: Model 2670 Controller Emission Indicator

Safety Interlock Keyswitch

The safety interlock keyswitch on the Model 2670 controller (Figure 2-4) limits laser access to only those personnel with permission and who have had adequate laser safety training. Without the key, the laser cannot be turned on. Nor can the laser be left on and the key removed. The key can only be removed in the "off" position. Always remove the key when done operating the laser.

Shutter

The laser head shutter is a lever-actuated mechanical device (Figure 2-3) that, when closed, interrupts the intracavity beam and prevents the laser from lasing. "O" is the open position, "Ø" is the closed position.

Cover Safety Interlocks

The Stabilite 2017 has safety interlocks for both the laser head cover and power supply cover. Removing either of these covers causes the main power contactor to open, which shuts off electrical power to the laser. The covers must be in place or their interlocks defeated before the laser will operate.

Danger!

Both the laser head and power supply contain electrical circuits operating at lethal voltage and current levels. Be extremely careful whenever the cover is removed from either unit. Avoid contact with the high voltage terminals and components inside each unit.

Laser Head

While the laser head cover is removed, be extremely careful to avoid exposure to laser or collateral radiation.

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